

- 6.1 The above concentration values and instrument conditions are for a Perkin-Elmer HGA- 2100, based on the use of a 20 μL injection, continuous flow purge gas and pyrolytic graphite.
- 6.2 Background correction may be required if the sample contains high dissolved solids.
- 6.3 The use of nitrogen as a purge gas is not recommended.
- 6.4 For every sample matrix analyzed, verification is necessary to determine that method of standard addition is not required (see part 5.2.1 of the Atomic Absorption Methods section of this manual).
- 6.5 If method of standard addition is required, follow the procedure given earlier in part 8.5 of the Atomic Absorption Methods section of this manual.

7.0 Precision and Accuracy

- 7.1 Precision and accuracy data are not available at this time.